



Thank you for choosing a Mediline product please read warranty and maintenance of your product.

Warranty applies from date of installation provided that installation is no more than 1 month from date of delivery. The warranty period is for 24 months against parts and labour, faulty workmanship and defective material on goods manufactured by Mediline.

Plus a further 12 months warranty on refrigeration system only.

Request for warranty are to be in writing and state model, date of installation, serial number and problem.

The warranty expressly excludes:

- Replacement of any glass or lights
- Any defects deemed by us to be attributed to negligence, misuse and/or accident in observing operating instructions
- Electrical faults and surges due to connection to inadequate power
- Loss of stock due to equipment malfunction **MEIDLINE WILL NOT ACCEPT ANY COSTS FOR STOCK LOSS** during warranty period under any circumstances.
- Electronic controller
- Compressor failure due to insufficient ventilation around equipment
- **Replacement Batteries.**



Set Up Instructions for Vaccine Refrigerators / Freezers

Dear Customer,

Thank you for choosing our refrigerator, to ensure that cabinet is installed correctly please follow the following steps.

- a. Unpack the refrigerator and place in position
- b. Ensure that refrigerator is level, if not you may adjust front levelling feet only
- c. Ensure there is 60mm ventilation space at rear and 30mm on sides of refrigerator for proper air flow.
- d. Position shelves at the desired levels that you require.
- e. Plug refrigerator into your 240 volts outlet – **TURN SWITCH ON AT REAR OF REFRIGERATOR** and let run for 2 hours during the period the refrigerator would have cycled a number of times. You are now ready to place your products on shelves.
- f. Cabinet temperature will vary form when the motor stops and starts. This is only air temperature and not product temperature.
- g. Temperature reading should only be done as follows, reset controller at end of day and do not open door again, next morning take your high and low read out before you open door.

Preventative Maintenance



It is important that condenser on back of cabinet is cleaned every 3 months, simply by brushing condenser in a vertical motion with a small brush or vacuum cleaner. However this will also depend on how dusty the area is. You may be able to do it every 6 months.


Refrigeration is fan forced to even ensure temperature throughout the cabinet with auto defrost and water disposal.



MEDI LINE[®]
A U S T R A L I A

Check Min/Max

Press  &  keys at same time for 3 seconds

000 will display – press  to go to 005 - Press √ once (controller unlocked)

Press √ for 3 seconds

$\bar{\Pi}$ Π II will display
(this is max temp)
Press √
Press √

Scroll up to find $\bar{\Pi}$ I n
(this is min temp)
Press √
Press √

How to Rest Min / Max

Scroll up to find CLr – Press √
Scroll to find 0 - Press √

Press √ for 3 seconds to go back to current temp.



MEDI LINE[®]
A U S T R A L I A

Alarm set points

Low alarm -50°C

High alarm -30°C

.

Delay on alarm is 10 minutes

Temperature set point


-40.0 °C



MEDI LINE[®]
A U S T R A L I A

HOW TO DOWNLOAD DATA USING USB

- Insert Data Logger into USB Slot on front of refrigerator

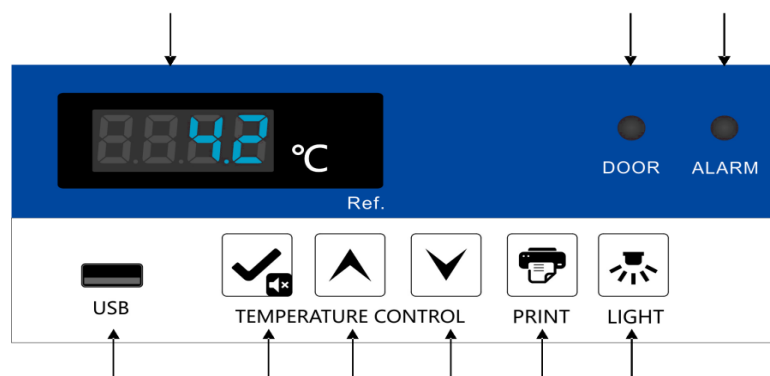

- Digital display will show **ON** to indicate data is downloading
- When finished display will show **END** – wait until display shows current temperature then remove USB
- Download information on to your computer and save as required.


Please note – Refrigerator will store the data for 2 months. At the start of each month the previous month will delete.

EG – On the 1st of March – Januarys data will delete and you will only be able to see February and March Data



MEDI LINE[®]
AUSTRALIA

How to change data logging recording time

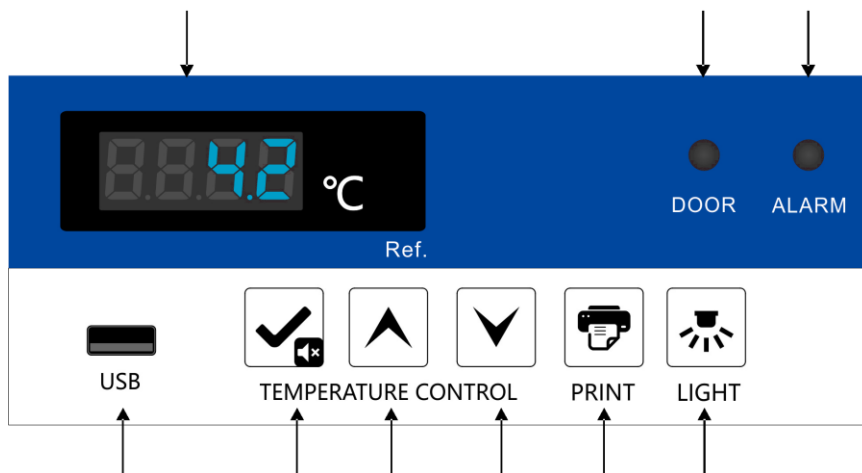

1. Press and hold Up & Down arrows at the same time until 000 is displayed
2. Use up arrow to go to 5, press the tick button
3. Hold the tick button until 000 is displayed
4. Use up arrow to go to 22, press the tick button
5. Press the down key until SCY
6. Press the tick button
Change to desired time EG 1 = 1 minute recording interval / 5 = 5-minute recording interval, then press the tick button
7. Press and hold tick button until refrigerators start beeping and display says LOF
8. Press and hold tick button and up arrow until refrigerator beeps and LOF is cleared from the screen

laftech
Contamination Control Solutions

Call: 1300 306 002 www.laftech.com.au